

<u>Title</u>	<u>Name</u>	<u>Approval Date</u>
Controls Infrastructure Group Leader	Ruslan Kadyrov	09/16/2015
ES&H Operations Manager	Lori Stiegler	09/16/2015
Quality Assurance Engineer	Joseph Zipper	09/16/2015

Serial No	Part No	Part Rev	ECN	Rev	ECN	Rev
	Beamline: <u>18ID - FXI</u>					
Deviation & Waiver: _____						

OP	Description	Name/Life #	Date	DR
10	Follow the ES&H and Personal Protective Equipment Requirements for the area.	H. Bassam 25077	10/10/17	

- 20 Verify measuring and test equipment used for this procedure contains a valid calibration label in accordance with NSLS-II Calibration Procedure PS-QAP-0901, where applicable.

H. Bassam 25077	10/10/17	
--------------------	----------	--

The technician is responsible for notifying the technical supervisor and/or the cognizant engineer of any discrepancies occurring during the performance of this procedure. All discrepancies shall be identified and reported in accordance with NSLS-II Discrepancy Reporting Procedure PS-QAP-0002.

- 30 BEAMLINE INFORMATION - This step shall be performed by the cognizant EPS Engineer.

G. Bischof 25835	10/10/17	
---------------------	----------	--

- A) Record the relevant Beamline name on this sheet, in the box for "Part No".
- B) Review this entire traveler and write in the relevant drawing, software, and procedure numbers where required

- 40 INSTALLATION VERIFICATION

- A) List the relevant drawing(s):

Drawing No.: LT-EL-BL-EI-EP-1180-90 Rev. No.: A

Drawing No.: _____ Rev. No.: _____

H. Bassam 25077	10/10/17	
--------------------	----------	--

- B) Verify the following items are acceptable:

- System layout configured per drawing(s)
- System Labeled / Tagged



OP	Description	Name/Life #	Date	DR
50	SOFTWARE CONFIGURATION A) Download PLC software to controller and verify it was accepted by the controller. B) Record the software part number: Part No.: <u>LT-EL-BL-EI- EPS-1180-70</u> Rev. No.: <u>A</u>	<u>G. BISCHOF</u> <u>25835</u>	<u>10/10/17</u>	
60	ACCEPTANCE TESTING A) List relevant testing procedure <u>PS-R-XFD-EPS-</u> Procedure No.: <u>CHK-001</u> Rev. No.: <u>1</u> B) Verify acceptable completion of test procedure C) Attach test report to this traveler	<u>G. Bischof</u> <u>25835</u>	<u>10/10/17</u>	
65	WATER LEAK DETECTION SYSTEM TEST Verify water leak detection system, as designed for this beamline, functions correctly.	<u>G. Bischof</u> <u>25835</u>	<u>10/10/17</u>	
70	Verify All Traveler Operations Complete	<u>H. Birsan</u> <u>25677</u>	<u>10/10/17</u>	
80	REVISION HISTORY (This step is informational and does not require signoff) Rev - Description - Date A First Release 8/27/2014 B OP#65 added 9/16/15 Joe DeLong removed as approver Ruslan Kadyrov added as approver			



Subject:	Beamline Equipment Protection System Test Checklist						
Number:	PS-R-XFD-EPS-CHK-001	Revision:	1	Effective:	8/19/14	Page:	1 of 8

Beam Line: 18-ID FXI

Test Date: 9/25/17

EPS Engineer: Garrett Bischof

BL Group Leader: Wah-Keat Lee

Beam Line master spreadsheet: Copy of LT-R-XFD-CO-DR-FXI-002

Pre-test setup:

Connect PPS interface test box at beam line EPS/PPS interface connector.

The Beam Line Master Spreadsheet contains a comprehensive list of all EPS related signals. As this test plan is executed note the results in the “test results” column of this spreadsheet.

Test Set 1: Vacuum

Vacuum Section:

Starting conditions: pressure at or below acceptable limits, GVs open.

Simulate pressure rise (toward poor vacuum) by disabling the vacuum gauge controller channel. Ensure EPICS PVs enter proper alarm states, photon shutter closes¹ and vacuum section is isolated. Two local gate valves and one (or more, if it is required to isolate upstream section, e.g. section branching) upstream gate valve close. Record results for the following:

EPICS PV Alarm

EPS fault

Photon Shutter closes

GVx closes

GVy close:

GVz closes

Attempt to open the front end shutter and gate valves through EPICS.

Photon shutter and gate valves cannot be opened through EPICS (1)

¹ if the intensity of the beam in the section cannot cause damage to its valves, poor vacuum doesn't have to cause shutter close

Brookhaven National Laboratory/ Photon Sciences Directorate

Subject: Beamline Equipment Protection System Test Checklist

Number: **PS-R-XFD-EPS-CHK-001** Revision: **1** Effective: **8/19/14** Page: **2 of 8**

Enable vacuum gauge controller and ensure EPICS alarms clear. Open gate valves and photon shutter through EPICS.

Photon shutter and gate valves can be opened through EPICS (2)

Gauge name:	EPS fault:				(1)	(2)
	Shutter:	d/s GV:	u/s GV:	D u/s:		
XF:18IDA-VA{DP:US-CCG:1}	Good	Good	Good	NA	Good	Good
XF:18IDA-VA{DP:US-TCG:1}	NA	NA	NA	NA	Good	Good
XF:18IDA-VA{DP:DS-CCG:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{DP:DS-TCG:1}	NA	NA	NA	NA	Good	Good
XF:18IDA-VA{Mir:CM-CCG:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{Mir:CM-TCG:1}	NA	NA	NA	NA	Good	Good
XF:18IDA-VA{TP:1-CCG:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{TP:1-TCG:1}	NA	NA	NA	NA	Good	Good
XF:18IDA-VA{WPFS:1-CCG:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{WPFS:1-TCG:1}	NA	NA	NA	NA	Good	Good
XF:18IDA-VA{PBSL:1-CCG:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{PBSL:1-TCG:1}	NA	NA	NA	NA	Good	Good
XF:18IDA VA{Mono:DCM-CCG:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{Mono:DCM-TCG:1}	NA	NA	NA	NA	Good	Good
XF:18IDA-VA{PMFS:1-CCG:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{PMFS:1-TCG:1}	NA	NA	NA	NA	Good	Good
XF:18IDA-VA{Mir:TM-CCG:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{Mir:TM-TCG:1}	NA	NA	NA	NA	Good	Good
XF:18IDA-VA{MFS:1-CCG:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{MFS:1-TCG:1}	NA	NA	NA	NA	Good	Good
XF:18IDA-VA{BT:1-CCG:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{BT:1-TCG:1}	NA	NA	NA	NA	Good	Good
XF:18IDA-VA{BT:1-CCG:2}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{BT:1-TCG:2}	NA	NA	NA	NA	Good	Good

Subject:	Beamline Equipment Protection System Test Checklist		
Number:	PS-R-XFD-EPS-CHK-001	Revision:	1
		Effective:	8/19/14
			Page: 3 of 8

Simulate pressure rise (toward poor vacuum) by disabling the vacuum pump controller channel. Ensure EPICS PVs enter proper alarm states, photon shutter closes² and vacuum section is isolated. Two local gate valves and one upstream (or more, if it is required to isolate upstream section, e.g. section branching) gate valve close. Record results for the following:

EPICS PV Alarm

EPS fault

Photon Shutter closes

GVx closes

GVy closes

GVz closes

Attempt to open the front end shutter and gate valves through EPICS.

Photon shutter and gate valves cannot be opened through EPICS (1)

Enable vacuum pump controller and ensure EPICS alarms clear. Open gate valves and photon shutter through EPICS.

Photon shutter and gate valves can be opened through EPICS (2)

Repeat for each vacuum section on beam line. For the sections adjacent to Front end and End Station, also simulate poor vacuum signal from remote systems.

EPS fault:	Front end:	End station:
Photon Shutter closes	Good	NA
GV2 closes	Good	NA
DP:DS GV closes	Good	NA
GVz closes	NA	NA
Shutter and GVs cannot be opened	Good	NA

Simulate a poor vacuum EPS signal to front end, GV2 and Shutter should close, and not be allowed to open.

EPS vacuum fault:	Front end:
Photon Shutter closes	Good
GV2 closes	Good
Shutter and GVs cannot be opened	Good

² if the intensity of the beam in the section cannot cause damage to its valves, poor vacuum doesn't have to cause shutter close

Brookhaven National Laboratory/ Photon Sciences Directorate

Subject: Beamline Equipment Protection System Test Checklist

Number: PS-R-XFD-EPS-CHK-001 Revision: 1 Effective: 8/19/14 Page: 4 of 8

Gauge name:	EPS fault:				-1	-2
	Shutter:	d/s GV:	u/s GV:	D u/s:		
XF:18IDA-VA{DP:US-IP:1}	Good	Good	Good	NA	Good	Good
XF:18IDA-VA{DP:DS-IP:1}	Good	Good	Good	NA	Good	Good
XF:18IDA-VA{Mir:CM-IP:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{TP:1-IP:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{WPFS:1-IP:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{PBSL:1-IP:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{Mono:DCM-IP:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{PMFS:1-IP:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{Mir:TM-IP:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{MFS:1-IP:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{BT:1-IP:1}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{BT:1-IP:2}	Good	Good	Good	Good	Good	Good
XF:18IDA-VA{BT:1-IP:3}	Good	Good	Good	Good	Good	Good

Test Set 2: Water flow

Water flow Section:

Record initial flow through section with all valves fully open.

Slowly close supply valve and record the LOW and LOW LOW PV Alarm levels.

If the channel is associated with an EPS action XFD-EPOS, also register the flow at which the EPS fault occurs. Ensure cable disconnection results in EPS action.

Repeat for each water circuit on beam line.

Sensor name:	Nominal flow:	LOW:	LOLO:	EPS:
XF:18IDA-OP{PPS:1}F:1-I	1.6	NA	NA	PPS only
XF:18IDA-OP{PPS:1}F:2-I	1.6	NA	NA	PPS only
XF:18IDA-OP{EPS:1}F-I	0.62	0.5	0.4	Closes FE shutter
XF:18IDA-OP{EPS:2}F-I	0.64	0.5	0.4	Closes FE shutter
XF:18IDA-OP{ }F:Main-I	2.73	NA	NA	NA
XF:18IDA-OP{Chl}F:Pmp1-I	2.0	NA	NA	NA
XF:18IDA-OP{Chl}F:Pmp2-I	2.1	NA	NA	NA
XF:18IDA-OP{Chl}F:Pmp3-I	2.0	NA	NA	NA

Brookhaven National Laboratory/ Photon Sciences Directorate

Subject:	Beamline Equipment Protection System Test Checklist						
Number:	PS-R-XFD-EPS-CHK-001	Revision:	1	Effective:	8/19/14	Page:	5 of 8

Test Set 3: Thermal

For each temperature transducer ensure the temperature measurement reports expected value.

Attach an appropriate adaptor and transducer simulator to the temperature input of the EPS system (remote IO chassis or Armor Block). Raise the temperature above each of the HI and HIHI PV alarm limits and ensure the alarm is reported. If the transducer is associated with an EPS interlock, continue to raise the temperature until the EPS trip level is exceeded. Ensure the appropriate EPS mitigation process occurs. Ensure channel cable disconnection results in EPS action.

Sensor name:	Temperature:	HI:	HIHI:	EPS:
XF:18IDA-OP{FM:1}T:T-I	28.3	150	200	Closes FE shutter
XF:18IDA-OP{FM:1}T:B-I	28.3	150	200	Closes FE shutter
XF:18IDA-OP{FM:1}T:I-I	28.3	150	200	Closes FE shutter
XF:18IDA-OP{FM:1}T:O-I	28.4	150	200	Closes FE shutter
XF:18IDA-OP{Mir:CM}T:Ax:X1-I	22.4	NA	NA	NA
XF:18IDA-OP{Mir:CM}T:Ax:X2-I	20.0	NA	NA	NA
XF:18IDA-OP{Mir:CM}T:1-I	20.0	70	80	Closes FE shutter
XF:18IDA-OP{Mir:CM}T:2-I	22.7	70	80	Closes FE shutter
XF:18IDA-OP{Mir:CM}T:CoolPlate1-I	20.1	70	80	Closes FE shutter
XF:18IDA-OP{Mir:CM}T:CoolPlate2-I	23.1	70	80	Closes FE shutter
XF:18IDA-OP{Mir:CM}T:HoldPlate1-I	22.3	70	80	Closes FE shutter
XF:18IDA-OP{Mir:CM}T:HoldPlate2-I	22.5	70	80	Closes FE shutter
XF:18IDA-OP{Mir:CM}T:DMsk1-I	22.8	70	80	Closes FE shutter
XF:18IDA-OP{Mir:CM}T:DMsk2-I	NA	NA	NA	NA
XF:18IDA-OP{Mir:CM}T:CmptnShld-I	22.4	70	80	Closes FE shutter
XF:18IDA-OP{WPFS:1}T:Ax:Y1-I	28.4	NA	NA	NA
XF:18IDA-OP{WSTPM:1}T:T-I	28.4	150	200	Closes FE shutter
XF:18IDA-OP{WSTPM:1}T:B-I	28.4	150	200	Closes FE shutter
XF:18IDA-OP{WSTPM:1}T:I-I	28.5	150	200	Closes FE shutter
XF:18IDA-OP{WSTPM:1}T:O-I	28.6	150	200	Closes FE shutter
XF:18IDA-OP{PBSL:1}T:T-I	28.4	NA	NA	NA
XF:18IDA-OP{PBSL:1}T:B-I	28.5	NA	NA	NA
XF:18IDA-OP{PBSL:1}T:I-I	28.3	NA	NA	NA
XF:18IDA-OP{PBSL:1}T:O-I	28.4	NA	NA	NA
XF:18IDA-OP{Mono:DCM}T:Ax:Y2-I	14.4	NA	NA	NA
XF:18IDA-OP{Mono:DCM}T:Ax:Th2-I	3.9	NA	NA	NA
XF:18IDA-OP{Mono:DCM}T:Ax:Chi2-I	-7.2	NA	NA	NA
XF:18IDA-OP{Mono:DCM}T:CmptnShld-I	-185.8	NA	NA	NA
XF:18IDA-OP{Mono:DCM}T:Crys1_1-I	-188.2	-150	-120	Closes FE shutter
XF:18IDA-OP{Mono:DCM}T:Crys1_2-I	-187.0	-150	-120	Closes FE shutter
XF:18IDA-OP{Mono:DCM}T:Crys2_1-I	-187.0	-150	-120	Closes FE shutter

Brookhaven National Laboratory/ Photon Sciences Directorate

Subject: Beamline Equipment Protection System Test Checklist

Number: PS-R-XFD-EPS-CHK-001 Revision: 1 Effective: 8/19/14 Page: 6 of 8

XF:18IDA-OP{Mono:DCM}T:Crys2_2-I	-185.4	-150	-120	Closes FE shutter
XF:18IDA-OP{Mono:DCM}T:Chi1Stage-I	17.1	NA	NA	NA
XF:18IDA-OP{Mono:DCM}T:Chi2Stage-I	13.5	NA	NA	NA
XF:18IDA-OP{PMFS:1}T:Ax:Y1-I	28.1	NA	NA	NA
XF:18IDA-OP{PBST:1}T:T-I	28.9	150	200	Closes FE shutter
XF:18IDA-OP{PBST:1}T:B-I	28.5	150	200	Closes FE shutter
XF:18IDA-OP{Mir:TM}T:Ax:X1-I	24.3	NA	NA	NA
XF:18IDA-OP{Mir:TM}T:Ax:X2-I	24.7	NA	NA	NA
XF:18IDA-OP{Mir:TM}T:HoldPlate1-I	24.7	NA	NA	NA
XF:18IDA-OP{Mir:TM}T:HoldPlate2-I	24.2	NA	NA	NA
XF:18IDA-OP{PCW}T:Sup-I	12.6	NA	NA	NA
XF:18IDA-OP{PCW}T:Rtn-I	13.8	NA	NA	NA
XF:18IDA-OP{DI:Water}T:Sup-I	28.7	NA	NA	NA
XF:18IDA-OP{DI:Water}T:Rtn-I	28.6	NA	NA	NA

Brookhaven National Laboratory/ Photon Sciences Directorate			
Subject:	Beamline Equipment Protection System Test Checklist		
Number:	PS-R-XFD-EPS-CHK-001	Revision: 1	Effective: 8/19/14 Page: 7 of 8

Test Set 4: EPS/PPS interface

For the primary Photon Shutter:

Front End: If **FRONT END ENABLE STATUS** and **FOE SECURE** and not **USER INTERLOCK** command the front end shutter to open.

Front end opens and **FRONT END OPEN STATUS** = OPEN:

Command the front end shutter to close.

Front end closes and **FRONT END OPEN STATUS** = CLOSED:

With the front end open force a user interlock.

Front end closes and **FRONT END OPEN STATUS** = CLOSED:

Attempt to open the front end shutter through EPICS while not all gate valves in white-beam region down to next secondary shutter are open.

Front end cannot be opened through EPICS:

For each secondary photon shutter:

If **PHOTON SHUTTER x ENABLE STATUS = 1 (ABILITY TO OPEN)** and **ESEE x SECURE** command the shutter to open.

Shutter opens and **PHOTON SHUTTER x OPEN STATUS** = OPEN

Command the shutter to close.

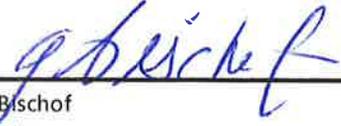
Shutter closes and **PHOTON SHUTTER x OPEN STATUS** = CLOSED

Attempt to open the shutter through EPICS while not all gate valves in white-beam region downstream the shutter are open.

Shutter cannot be opened through EPICS:

Brookhaven National Laboratory/ Photon Sciences Directorate			
Subject:	Beamline Equipment Protection System Test Checklist		
Number:	PS-R-XFD-EPS-CHK-001	Revision: 1	Effective: 8/19/14 Page: 8 of 8

Prepared by:

X 

Garrett Bischof
Document Preparer

Approved by:

X 

Harmanpreet Bassan
Deputy Controls Group Leader

X 

Wah-Keat Lee
Beamline Group Leader